Levenhuk

Digital

EN	User M
CZ	Návod
DE	Bedie
ES	Guía d
PL	Instru
RU	Инстр

230x
ZOOM

2

400x
ZOOM

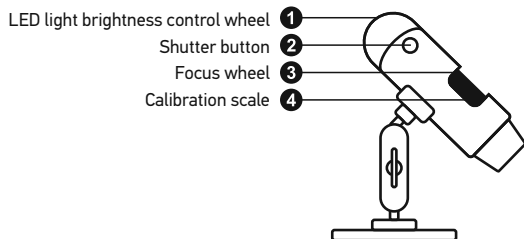
1.3

leve

Levenhuk DTX 30 DTX 50

Digital Microscope

EN

The kit includes: microscope, Levenhuk DTX installation CD, calibration scale, user guide.

Using the microscope

Remove the protective cap from the objective lens. Use the focus wheel to focus your view. The shutter button allows you to capture images with default settings (using MicroCapture software to create images with default settings is described in the corresponding section of this User Guide).

If you wish to adjust brightness of the LED illumination, you can use the illumination brightness adjustment ring at the top of the microscope body.

Make sure the microscope is connected to your PC before proceeding with the setup. Insert the installation CD into your CD drive. The Main Menu window will be displayed. Click the **install Microscope Driver** to install the required drivers and follow the instructions on screen. After the drivers are installed, click the **MicroCapture software** to install the software. Keep in mind that a system restart will be required during installation. The installation CD contains a copy of this User Guide, which you can read by clicking the corresponding entry in the Main Menu.

Never, under any circumstances, look directly at the sun or another bright source of light or as this may cause **PERMANENT** eye damage and may lead to **BLINDNESS**.

Image resolution

From the **Options** menu, select **Preview Size**.

Rotating the image

From the **Options** menu, select **Rotate** and choose clockwise or counterclockwise.

Date and time

From the **Options** menu, select **Date/Time** to set the date and time during preview.

Language

From the **Options** menu, select **Language** to choose the language.

Full screen mode

There are three ways to enter full screen mode:

1. Double-click an image in the main window;
2. Click the **Full screen mode** icon on the toolbar;
3. From the **Options** menu, select **Fullscreen**.

To exit full screen mode, click the corresponding icon on the toolbar.

Skins

From the **Options** menu, select **Skin** to change the skin.

Settings

From the **Options** menu, select **Preferences** to set the application, including brightness, contrast and gamma.

Capturing an image

Click the corresponding icon on the toolbar or press F11 on your keyboard. A thumbnail of the captured image will appear on the side of the main window.

Recording a video

Click the corresponding icon on the toolbar or select **Video** from the **Capture** menu. The camcorder icon on the toolbar will change once you start recording. To stop recording you can either click the corresponding icon on the toolbar again or select **Stop capturing** from the **Capture** menu.

Note: recommended video resolution is 640x480.

Preview

Left-click the thumbnail of the image you wish to edit and it will open in the main window of the application. The image resolution will be displayed in the bottom left corner of the screen. Double-clicking a thumbnail will open the image in a separate window.

Saving images

To save an image, right-click the thumbnail of the desired image and select **Save** from the drop-down menu. Images may be saved in .jpeg and .bmp formats. When saving images in .jpeg format, clicking **Advanced** allows you to view and edit additional options.

Deleting images

To delete an image, right-click the thumbnail of the desired image and select **Delete** from the drop-down menu. When viewing images in a separate window, you can click the corresponding icon on the toolbar to delete the image.

Copying images

To copy an image, right-click the thumbnail of the desired image and select **Copy** from the drop-down menu.

Calibration and measurements

Place a ruler on the stage, focus your view and aimage. Take note of the current magnification by looking at the calibration scale on the microscope body. Double-click the image thumbnail to open it in a separate window. Enter the magnification value into the corresponding field in the upper right corner of the preview window. Upon calibrating the microscope, you will be able to measure various parameters of observed specimens. All the buttons on the toolbar of the preview window are described below.

Line

Press and hold the left mouse button to start drawing a line. Release the left mouse button to complete the line.

Free-form line

Press and hold the left mouse button and draw a line. Release the left mouse button to complete the segment. You can continue adding more segments to the free-form line.

Radius

Draw a straight line on the image. The application will calculate the radius, as well as circumference and area of the corresponding circle.

Diameter

Draw a straight line on the image. The application will calculate the diameter, as well as circumference and area of the corresponding circle.

Angle

Press and hold the left mouse button to draw a line. Release the left mouse button and draw another line. The application will calculate the angle between the two lines.

Measurement units

From the drop-down menu on the toolbar you can select the units for calculations (pixels, inches, kilometers, meters, etc.).

Press this button to create a line.

Press this button to edit the line.

These buttons allow you to draw a line.

Press this button to choose a color.

Press this button to choose a line style.

Press this button to choose a line width.

Press this button to undo the last action.

Note:

1. Magnification range of Levenhuk DTX 50 is 200x-1000x. Do not use magnification values that exceed this range.
2. The application calculates values based on the image resolution. The results may not be precise.
3. Measurement functions are only available to the user with the appropriate permissions.

Working with the application for Mac OS users

Installing software

Browse to the **mac** folder on the installation CD and copy **MicroCapture** file to the desktop.

Operating the microscope

Use a USB cable to connect the microscope to your computer. Run the MicroCapture application by double-clicking its icon.

Observing specimens

A magnified image of the observed specimen may be seen in the main window of the application.

Image resolution

From the **Options** menu, select **Preview Size** to set the image resolution. Keep in mind that the preview window cannot be resized (standard resolution is 640x480). The image resolution is shown in the bottom left corner of the preview window.

Date and time

From the **Options** menu, select **Date/Time** to display or hide the date and time of capture during preview.

Capturing an image

Click the corresponding icon on the toolbar or select **Photo** from the **Capture** menu to capture an image. An image thumbnail will appear on the right side of the main window.

Recording a video

Click the corresponding icon on the toolbar or select **Video** from the **Capture** menu to start recording a video.

Note: during recording, the camcorder icon on the toolbar will turn red. Click the icon again to stop recording.

Editing images and videos

Image and video thumbnails are displayed on the right side of the main window of the application. Double-click an image thumbnail to open it in a separate preview window. Editing images with MicroCapture application is similar to working with other applications in Mac OS.

Care and maintenance

Never, under any circumstances, look directly at the laser through this device, as this may cause PERMANENT EYE DAMAGE. BLINDNESS. White LEDs are very bright; do not look directly at them. Do not damage to your vision. Take necessary precautions. Do not use the device if you are others who have not read or who do not fully understand the instructions. Do not use your microscope and before using it for the first time, read the instructions for every component and connection. Do not try to repair or clean the device on your own. For repairs and cleaning of any kind, please contact the manufacturer. Do not center. Protect the device from sudden impact or vibration. Do not apply excessive pressure when adjusting focus. Do not touch the device. Do not touch the optical surfaces with your fingers. To clean the device, use only soft, lint-free wipes and special optics cleaning tools from Leica. Do not use the device for lengthy periods of time, and do not leave the device away from water and high humidity. Do not use the device to replace the dust cover after you are finished with the device. Do not use the device and stains. If you are not using your microscope, store the objective lenses and eyepieces separately from the device. Do not place the device away from hazardous acids and other chemicals. Do not place the device near other sources of high temperatures. To avoid damage to the device, use the application first, and then disconnect the microscope from the power source. Do not illuminate goes off completely, and disconnect the device immediately if a small part or a battery is swollen or damaged.

Battery safety instructions

Always purchase the correct size and grade of batteries. Do not use old or worn batteries. Always replace the whole set of batteries at once, or batteries of different types. Clean the device before battery installation. Make sure the battery is inserted with the correct polarity (+ and -). Remove batteries from equipment after use. Do not use equipment for a long period of time. Remove used batteries promptly after use, as this may cause leakage, fire, or explosion. Do not use the device at high temperatures, leakage, or explosion. Never use the device near fire. Remember to switch off devices after use. Keep the device away from children. Do not risk of ingestion, suffocation, or poisoning.

Specifications

Image sensor	1.3Mpx, interpolated to 2Mpx (DTX 50) 2Mpx, interpolated to 5Mpx (DTX 30)
Magnification	20x–400x (DTX 50), 20x–230x (DTX 30)
Available image resolutions (DTX 50)	1600x1200, 1280x1024, 1280x960, 1024x768, 800x600, 640x480, 352x288, 320x240, 160x120
Available image resolutions (DTX 30)	2560x2048 (5M), 2000x1600, 1600x1280 (2M), 1280x1024, 1024x960, 1024x768, 800x600, 640x480, 352x288, 320x240, 160x120
Available video resolutions (DTX 50)	1600x1200, 1280x1024, 1280x960, 1024x768, 800x600, 640x480, 352x288, 320x240, 160x120
Available video resolutions (DTX 30)	2560x2048 (5M), 2000x1600, 1600x1280 (2M), 1280x1024, 1024x960, 1024x768, 800x600, 640x480, 352x288, 320x240, 160x120
Photo/video format	*.jpeg or *.bmp/*.avi
Focusing	Manual, 0–150mm focusing range
Illumination	8-LED system with variable brightness
Power source	USB 2.0 cable
Frame rate	30fps
Software languages	English, German, Spanish, Korean, French, Russian
Dimensions (length x diameter)	125x33mm

Levenhuk reserves the right to modify or discontinue any product without prior notice.

System requirements

Operating system: Windows XP SP2/Vista/7/8, Mac 10.6–10.8,
CPU: at least P4 1.8GHz, RAM: 512MB, interface: USB 2.0

Levenhuk International Lifetime Warranty

All Levenhuk telescopes, microscopes, binoculars and other optical products, except for accessories, carry a lifetime warranty against defects in materials and workmanship. Lifetime warranty is a guarantee on the lifetime of the product on the market. Levenhuk accessories are warranted to be free of defects in materials and workmanship for six months from date of retail purchase. Levenhuk will repair or replace such product or part thereof which, upon inspection by Levenhuk, is found to be defective in materials or workmanship. As a condition to the obligation of Levenhuk to repair or replace such product, the product must be returned to Levenhuk together with proof of purchase satisfactory to Levenhuk. This warranty does not cover consumable parts, such as bulbs (electrical, LED, halogen, energy-saving and other types of lamps), batteries (rechargeable and non-rechargeable), electrical consumables etc.

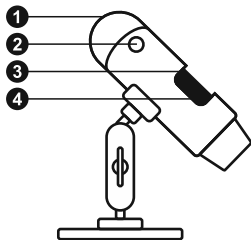
For further details, please visit our web site: www.levenhuk.com/warranty

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Levenhuk® is a registered trademark of Levenhuk, Inc. © 2006-2015 Levenhuk, Inc. All rights reserved.
Levenhuk, Inc. 924-D East 124th Ave. Tampa, FL 33612 USA